
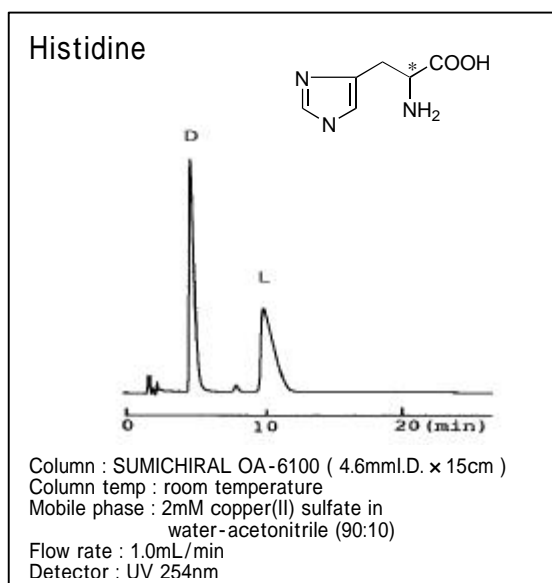
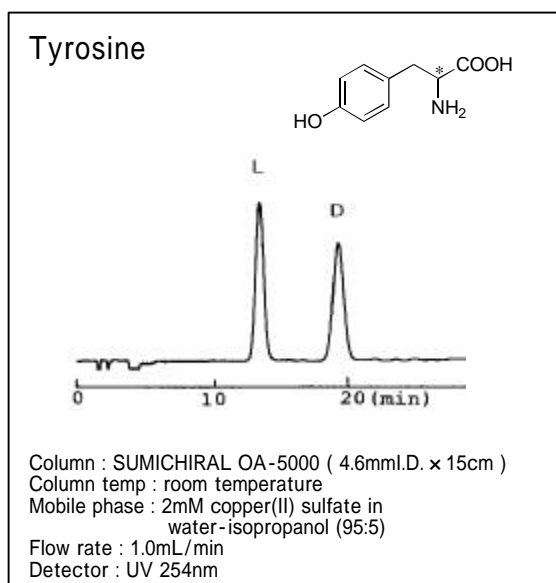
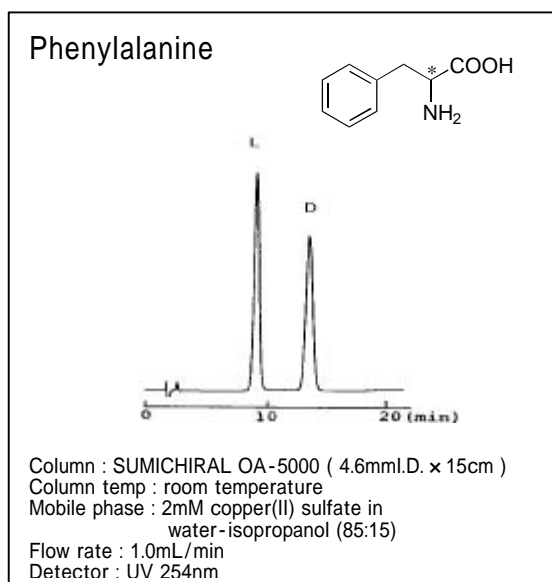


SUMICHIRAL TECHNICAL NEWS

(Vol. 9 7 0 9)

Enantiomer Separation of Aromatic Amino Acids by SUMICHIRAL OA

SUMICHIRAL OA-5000 and OA-6100 are chiral ligand exchange HPLC columns, and are extremely effective for direct enantiomer separation of wide range of amino acids. In this report, we introduce our applications of enantiomer separation of aromatic amino acids.

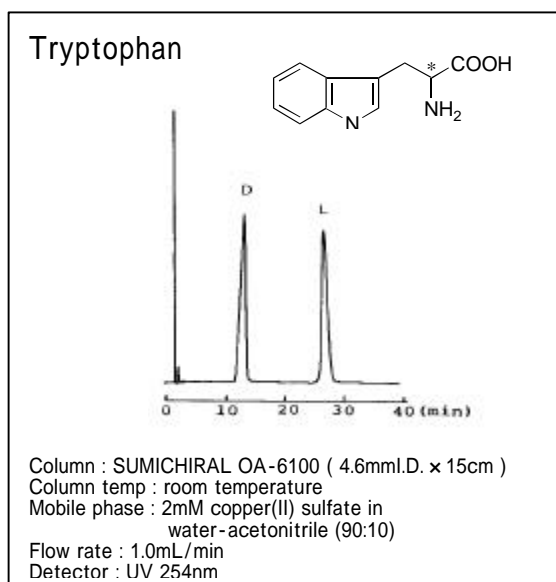

SUMIKA CHEMICAL ANALYSIS SERVICE, Ltd.
Scientific Instruments Div. Chromatography Group

