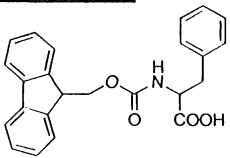
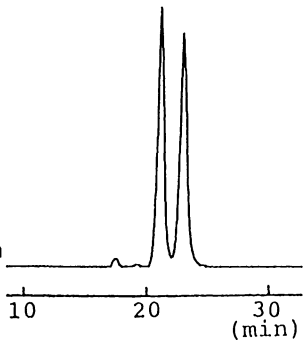
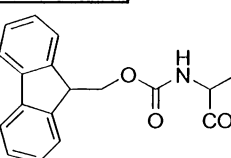
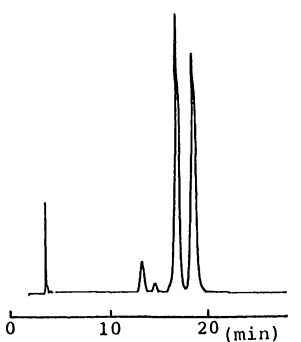
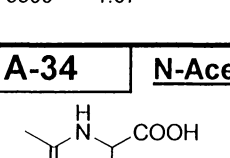
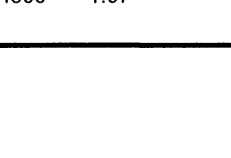
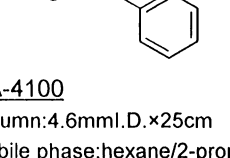
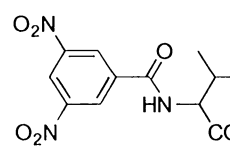
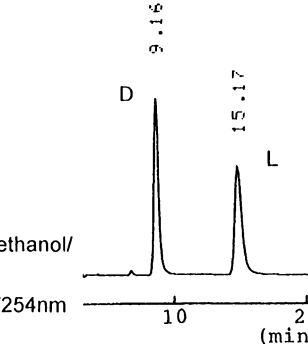
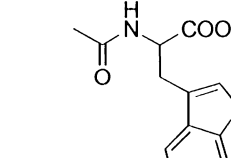
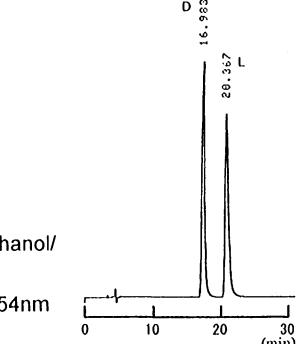


A-30  <u>OA-3200</u> Column:4.6mmI.D.×25cm Mobile phase:0.01mol/L ammonium acetate in methanol Flow rate:1.0mL/min Detector:UV254nm	N-FMOC-phenylalanine 	A-31  <u>OA-3200</u> Column:4.6mmI.D.×25cm Mobile phase:0.01mol/L ammonium acetate in methanol Flow rate:1.0mL/min Detector:UV254nm	N-FMOC-valine 
OA-3200 1.10 OA-3300 1.08 OA-7000 1.05		OA-3200 1.12 OA-3300 1.08	
A-32  <u>OA-3200</u> Column:4.6mmI.D.×25cm Mobile phase:0.01mol/L ammonium acetate in methanol Flow rate:1.0mL/min Detector:UV254nm	N-FMOC-leucine 	A-33  <u>OA-4500</u> Column:4.6mmI.D.×25cm Mobile phase:hexane/2-propanol/methanol/trifluoroacetic acid (90:5:5:0.2) Flow rate:1.0mL/min Detector:UV254nm	N-Formylphenylalanine
OA-3200 1.20 OA-3300 1.07		OA-4700 1.12 OA-4500 1.07	
A-34 N-Acetylphenylalanine  <u>OA-4100</u> Column:4.6mmI.D.×25cm Mobile phase:hexane/2-propanol/methanol/trifluoroacetic acid (90:5:5:0.2) Flow rate:1.0mL/min Detector:UV254nm Injection:5μL (5mg/mL)			
OA-4700 1.38 { D, L } OA-4700 1.24 { D, L } OA-4100 1.18	OA-4600 1.15 OA-4600 1.09 { D, L } OA-4000 1.08	OA-4500 1.08	
A-35  <u>OA-4800</u> Column:4.6mmI.D.×25cm Mobile phase:hexane/2-propanol/methanol/trifluoroacetic acid (70:20:10:0.2) Flow rate:1.0mL/min Detector:UV254nm Injection:0.5μL (2mg/mL)	N-3,5-Dinitrobenzoylvaline 	A-36  <u>OA-4700</u> Column:4.6mmI.D.×25cm Mobile phase:hexane/2-propanol/methanol/trifluoroacetic acid (90:5:5:0.2) Flow rate:1.0mL/min Detector:UV254nm Injection:1μL (10mg/mL)	N-Acetyltryptophan 
OA-4800 2.02 { D, L } OA-4900 1.30 { L, D }		OA-4700 1.25 { D, L }	