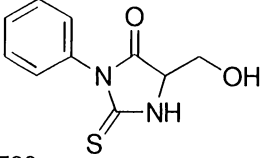
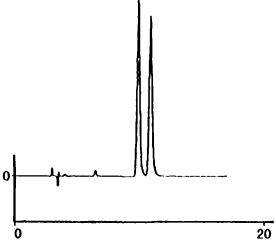
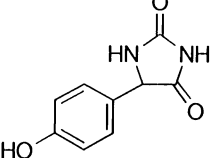
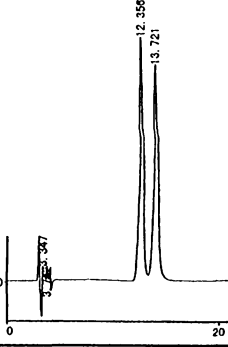
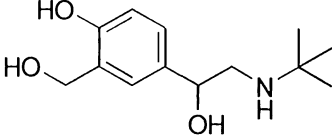
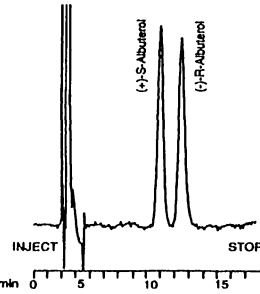
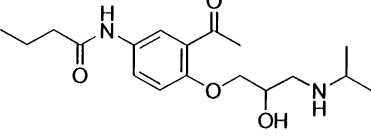
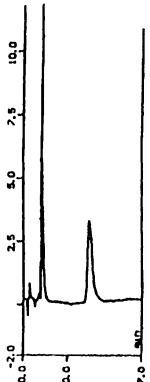
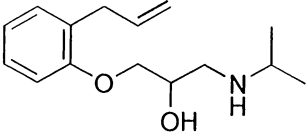
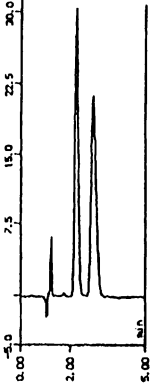
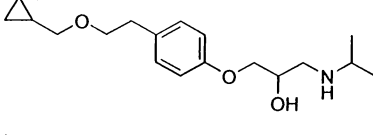
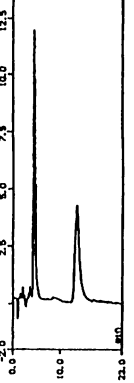
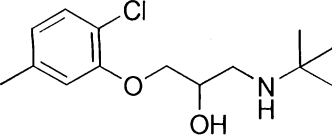
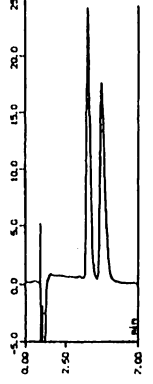
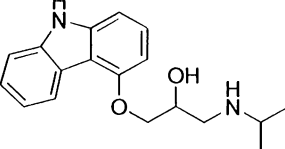
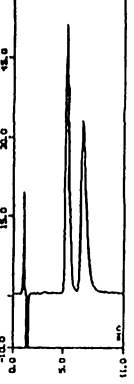


B-33 PTH-Serine  <u>OA-4700</u> Column: 4.6mm I.D. × 25cm Mobile phase: hexane/2-propanol/methanol/trifluoroacetic acid (80:15:5:0.2) Flow rate: 1.0mL/min Detector: UV254nm Injection: 1μL (1mg/mL)  OA-4700 1.23 OA-4100 1.06	B-34 5-(4-Hydroxyphenyl) hydantoin  <u>OA-4700</u> Column: 4.6mm I.D. × 25cm Mobile phase: hexane/2-propanol/methanol/trifluoroacetic acid (80:15:5:0.2) Flow rate: 1.0mL/min Detector: UV254nm Injection: 5μL (1mg/mL)  OA-4700 1.15
B-35 Albuterol  <u>OA-4900</u> Column: 4.6mm I.D. × 25cm Mobile phase: hexane/1,2-dichloroethane/methanol/trifluoroacetic acid (160:93:16:0.67) Flow rate: 1.0mL/min Detector: UV276nm  OA-4900 1.18 { +, - }	B-36 Acebutolol  <u>CBH</u> Column: 4.0mm I.D. × 10cm Mobile phase: 5% 2-propanol in 10mmol/L sod.ac.b. pH5.5 + 50μmol/L EDTA Flow rate: 0.9mL/min Detector: UV225nm  CBH 4.25
B-37 Alprenolol  <u>AGP</u> Column: 4.0mm I.D. × 10cm Mobile phase: 3% acetonitrile in 10mmol/L sod.ac.b. pH4.0 (total acetate conc. = 60mmol/L) Flow rate: 0.9mL/min Detector: UV225nm  AGP 1.61	B-38 Betaxolol  <u>CBH</u> Column: 4.0mm I.D. × 10cm Mobile phase: 5% 2-propanol in 10mmol/L sod.ac.b. pH5.5 + 50μmol/L EDTA Flow rate: 0.9mL/min Detector: UV225nm  CBH 3.37
B-39 Bupranolol  <u>AGP</u> Column: 4.0mm I.D. × 10cm Mobile phase: 0.5% 2-propanol in 40mmol/L amm.ac.b. pH4.1 (total acetate conc. = 190mmol/L) Flow rate: 0.9mL/min Detector: UV225nm  AGP 1.37	B-40 Carazolol  <u>AGP</u> Column: 4.0mm I.D. × 10cm Mobile phase: 0.5% 2-propanol in 4.9mmol/L amm.ac.b. pH4.1 (total acetate conc. = 25mmol/L) Flow rate: 0.9mL/min Detector: UV225nm  AGP 1.38