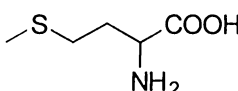
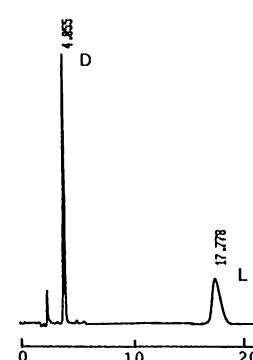
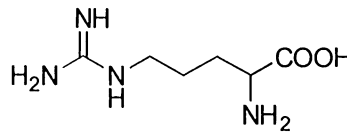
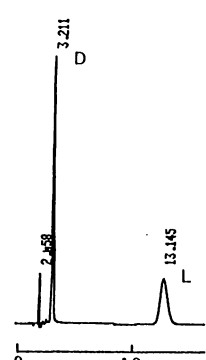
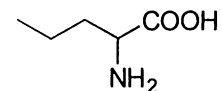
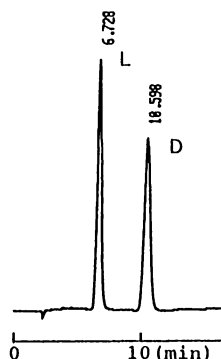
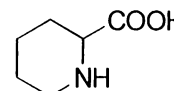
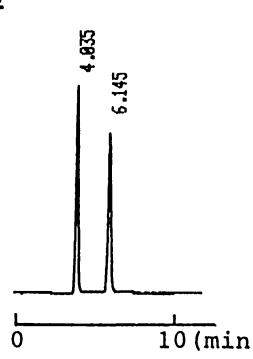
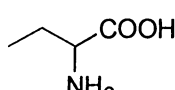
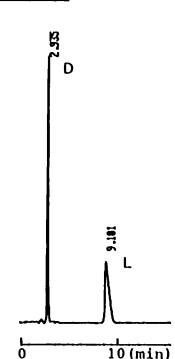
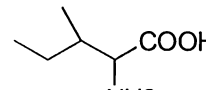
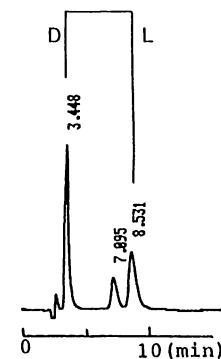
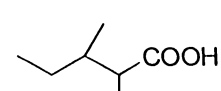
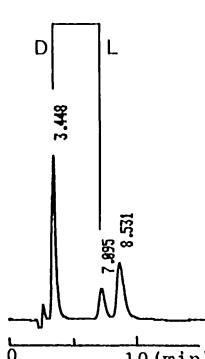
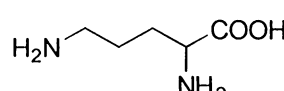
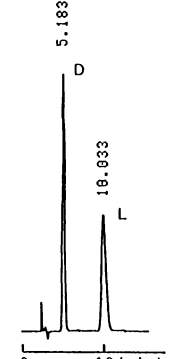


G-09 Methionine  <u>OA-6100</u> Column: 4.6mm I.D. × 15cm Mobile phase: 2mmol/L copper(II) sulfate in water/acetonitrile (95:5) Flow rate: 1.0mL/min Detector: UV254nm Injection: 5μL (1.3mg/mL)  OA-6100 6.16 { D, L } OA-6000 4.95 { D, L } OA-5500 1.30 { D, L } OA-5000 1.28 { L, D }	G-10 Arginine  <u>OA-6100</u> Column: 4.6mm I.D. × 15cm Mobile phase: 2mmol/L copper(II) sulfate in water/acetonitrile (98:2) Flow rate: 1.0mL/min Detector: UV254nm Injection: 1μL (5mg/mL)  OA-6100 6.50 { D, L } OA-6000 2.39 { D, L } OA-5000 2.15 { L, D }
G-11 Norvaline  <u>OA-5000</u> Column: 4.6mm I.D. × 15cm Mobile phase: 2mmol/L copper(II) sulfate in water/2-propanol (95:5) Flow rate: 1.0mL/min Detector: UV254nm Injection: 3μL (2mg/mL)  OA-6100 8.46 { D, L } OA-6000 5.44 { D, L } OA-5000 1.72 { L, D }	G-12 Pipecolic acid  <u>OA-5500</u> Column: 4.6mm I.D. × 15cm Mobile phase: 1mmol/L copper(II) sulfate in water Flow rate: 1.0mL/min Detector: UV254nm  OA-5500 1.80 OA-5000 1.77 { L, D } OA-6000 1.41 { L, D } OA-6100 1.14
G-13 2-Amino-n-butyric acid  <u>OA-6100</u> Column: 4.6mm I.D. × 15cm Mobile phase: 2mmol/L copper(II) sulfate in water/acetonitrile (98:2) Flow rate: 1.0mL/min Detector: UV254nm Injection: 1μL (5mg/mL)  OA-6100 5.00 { D, L } OA-6000 2.88 { D, L } OA-5000 1.80 { L, D }	G-14 Isoleucine  <u>OA-6000</u> Column: 4.6mm I.D. × 15cm Mobile phase: 2mmol/L copper(II) sulfate in water/acetonitrile (85:15) Flow rate: 1.0mL/min Detector: UV254nm  OA-6000 3.49 { D, L } OA-5000 1.72 { L, D } OA-6100 1.45 { D, L } OA-5500 1.15 { L, D }
G-15 Alloisoleucine  <u>OA-6000</u> Column: 4.6mm I.D. × 15cm Mobile phase: 2mmol/L copper(II) sulfate in water/acetonitrile (85:15) Flow rate: 1.0mL/min Detector: UV254nm  OA-6000 2.79 { D, L } OA-5000 1.67 { L, D } OA-5500 1.20 { L, D }	G-16 Ornithine  <u>OA-6100</u> Column: 4.6mm I.D. × 15cm Mobile phase: 1mmol/L copper(II) sulfate in water Flow rate: 1.0mL/min Detector: UV254nm Injection: 2μL (5mg/mL)  OA-6100 2.29 { D, L } OA-5000 1.38 { L, D }