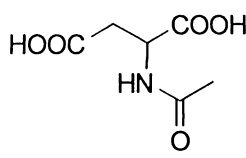
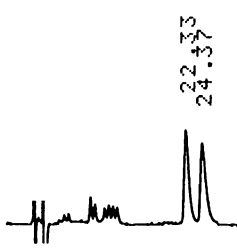
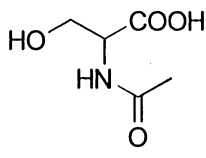
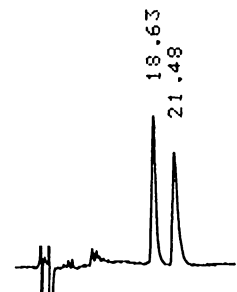
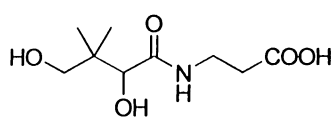
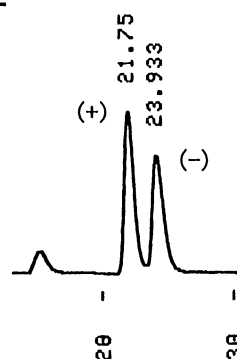
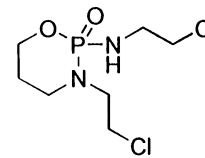
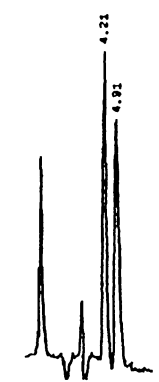
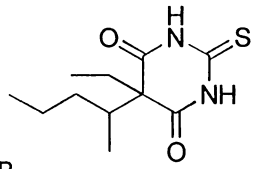
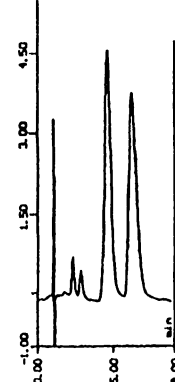
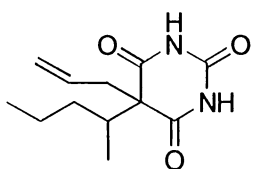
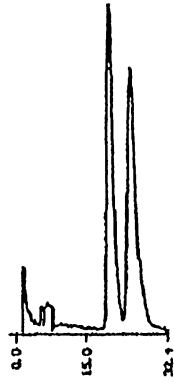
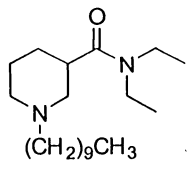
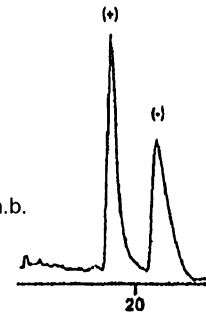
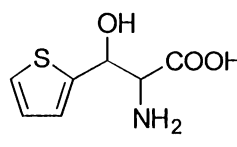
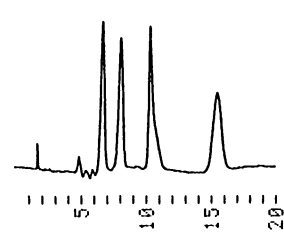


G-49 N-Acetyl-aspartic acid  <u>OA-4700</u> Column:4.6mmI.D.×25cm Mobile phase:hexane/2-propanol/methanol/ trifluoroacetic acid (90:5:5:0.2) Flow rate:1.0mL/min Detector:UV230nm Injection:2μL (10mg/mL)	 OA-4700 1.11	G-50 N-Acetylserine  <u>OA-4700</u> Column:4.6mmI.D.×25cm Mobile phase:hexane/2-propanol/methanol/ trifluoroacetic acid (90:5:5:0.2) Flow rate:1.0mL/min Detector:UV230nm Injection:2μL (10mg/mL)	 OA-4700 1.18
G-51 Pantothenic acid  <u>OA-6100</u> Column:4.6mmI.D.×15cm×2 Mobile phase:2mmol/L copper(II) sulfate in water/acetonitrile (95:5) Flow rate:1.0mL/min Detector:UV254nm	 OA-6100 1.12 { +, - }	G-52 Ifosfamide  <u>AGP</u> Column:4.0mmI.D.×10cm Mobile phase: 1% 2-propanol in 10mmol/L sod.ph.b. pH7.0 Flow rate:0.9mL/min Detector:UV195nm	 AGP 1.23
G-53 Penthiobarbital  <u>AGP</u> Column:4.0mmI.D.×10cm Mobile phase: 6% 2-propanol in 100mmol/L sod.ph.b. pH6.0 Flow rate: Detector:	 AGP 1.48	G-55 Secobarbital  <u>AGP</u> Column:4.0mmI.D.×10cm Mobile phase: 0.3% 2-propanol in 100mmol/L sod.ph.b. pH6.0 Flow rate:0.9mL/min Detector:UV225nm	 AGP 1.24
G-56 1-Decyl-3-(N,N-diethyl-carbamoyl)piperidine  <u>AGP</u> Column:4.0mmI.D.×10cm Mobile phase:10% ethanol in 25mmol/L sod.ph.b. pH7.0 + 25mmol/L tetrabutylammonium HSO4 Flow rate:0.9mL/min Detector:UV225nm	 AGP 1.56 { +, - }	G-57 3-(2-Thienyl)serine  <u>OA-5000</u> Column:4.6mmI.D.×15cm Mobile phase:2mmol/L copper(II) sulfate in water/methanol (70:30) Flow rate:1.0mL/min Detector:UV254nm	 OA-5000 OA-4800 OA-4900